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晉色園生物科技流動實驗室計劃
Sik Sik Yuen Biotechnology MobileLab Program

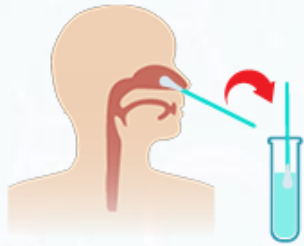
晉色園主辦可譽中學暨可譽小學
Ho Yu College and Primary School (Sponsored by Sik Sik Yuen)

Comparison of COVID-19 Molecular Tests using Nasopharyngeal (NP) specimen collection and SalivaDirect collection

Molecular Tests (Nucleic Acid Detection)

1. Obtain Specimen:
NP swab

2. Extract RNA
from specimen
and convert to
DNA.

3. Amplify by PCR
with SARS-CoV-2
specific primers.

4. Interpret results:
presence of viral
RNA indicates
active SARS-CoV-2
infection.

SalivaDirect Test

1. Obtain Specimen:
SalivaDirect

2. RNA-extraction free
process. Convert RNA from
specimen to DNA and
amplify by PCR with SARS-
CoV-2 specific primers.

3. Interpret results:
presence of viral
RNA indicates
active SARS-CoV-2
infection.

mRNA?
DNA?
PCR?

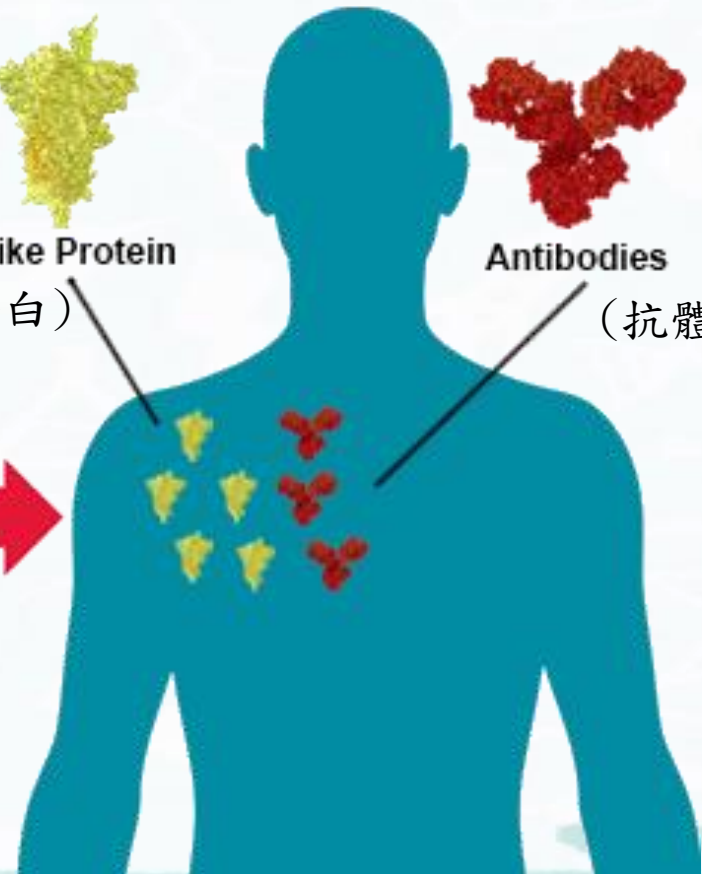
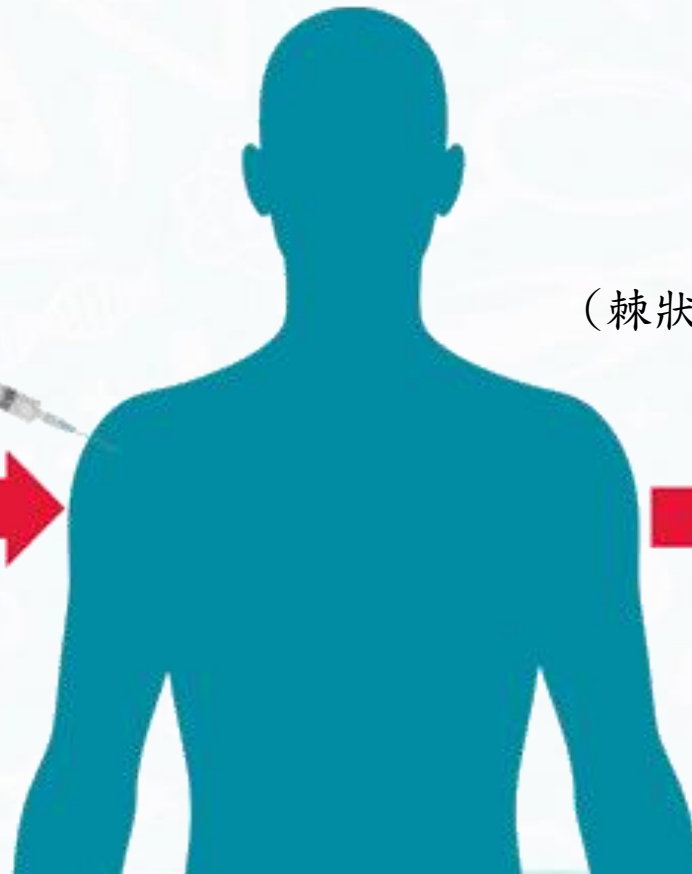
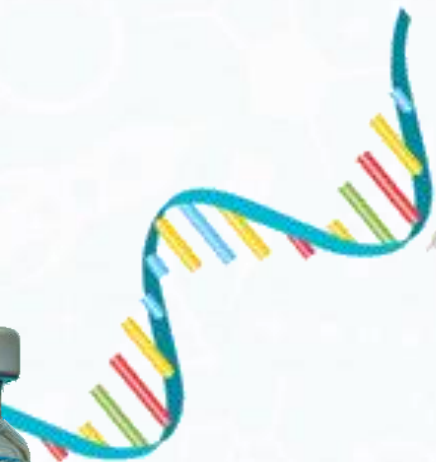
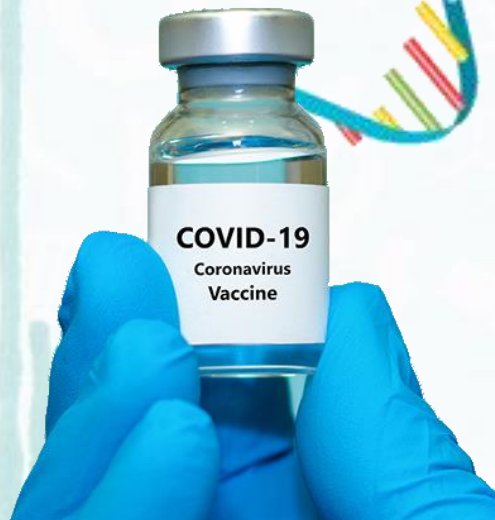
mRNA for spike protein of coronavirus

mRNA is injected into muscle


The vaccine triggers production of the spike protein and antibodies specific to it

Spike Protein
(棘狀蛋白)


Antibodies
(抗體)

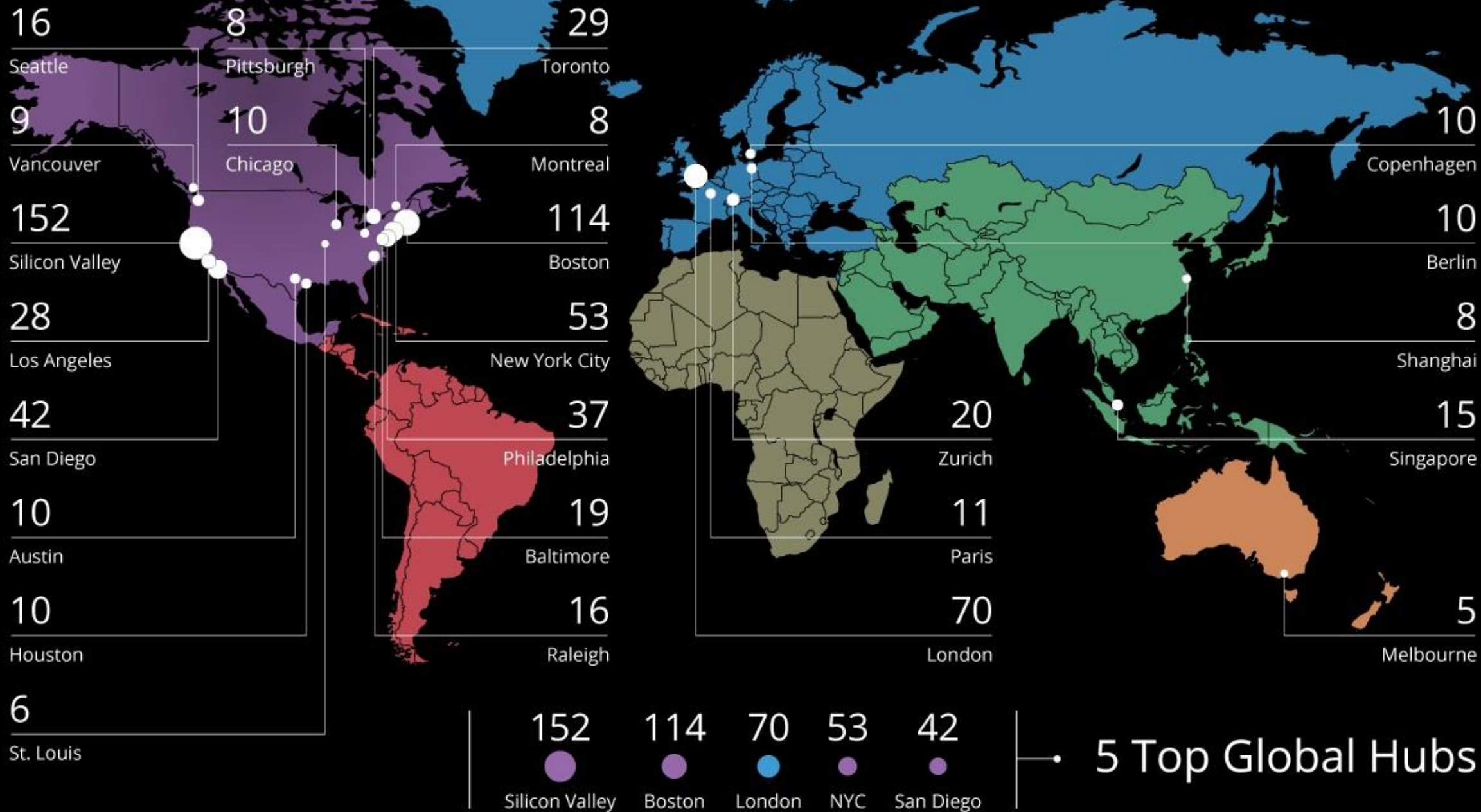Biotechnology


U.S. Biotechnology Market

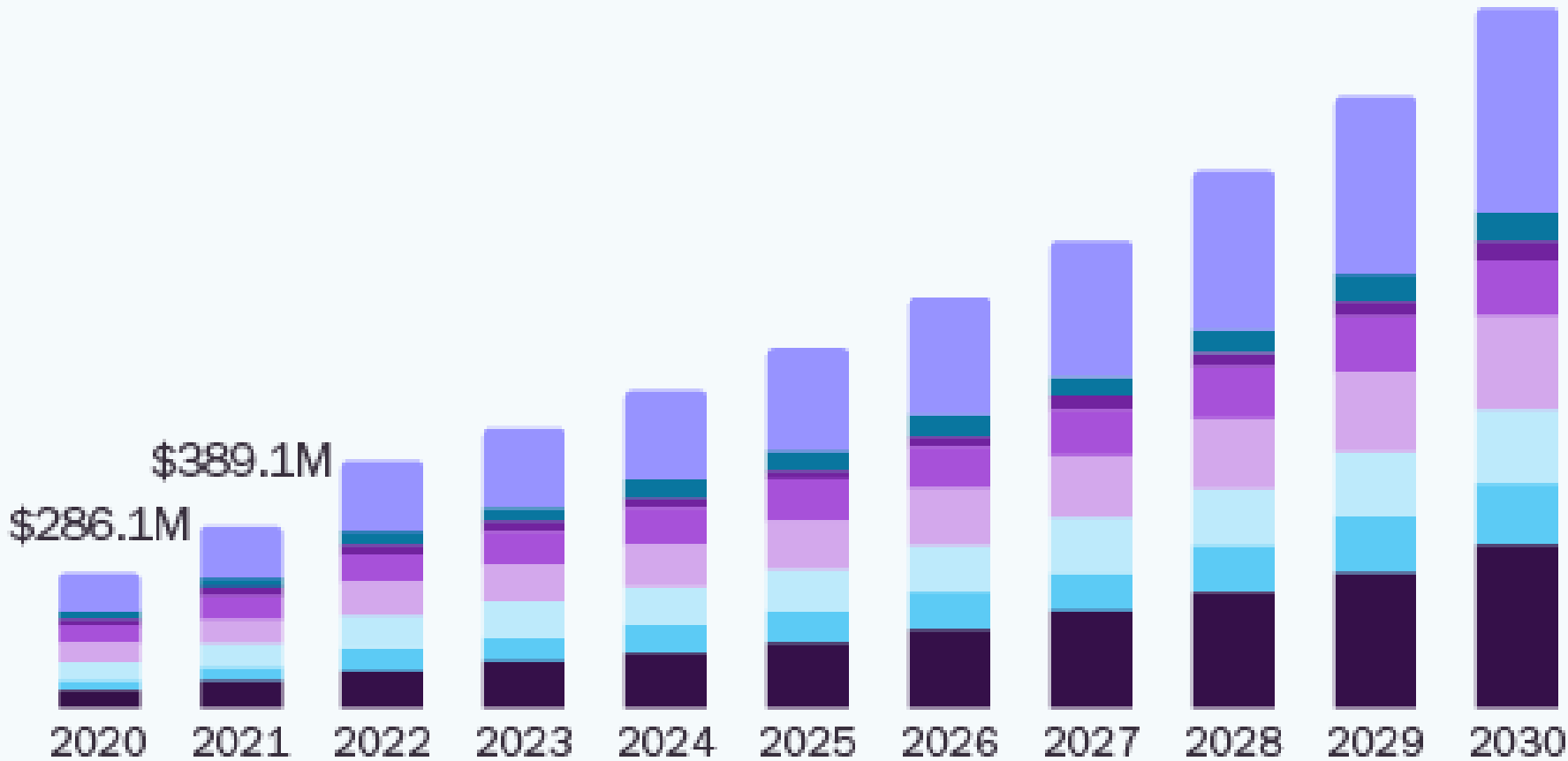size, by technology, 2020 - 2030 (USD Million)

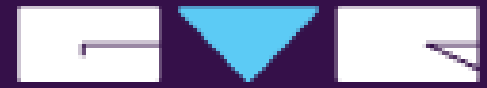- DNA Sequencing
- Nanobiotechnology
- Tissue Engineering & Regeneration
- Fermentation
- Cell-based Assays
- PCR Technology
- Chromatography
- Others

GRAND VIEW RESEARCH

13.9%

U.S. Market CAGR,
2022 - 2028

Source:
www.grandviewresearch.com

**Hands-on Biotech: The Lawrence Hall of Science
(University of California, Berkeley)**

Science Adventure Lab (2009 Launched)

North Carolina's BioNetwork Mobile Laboratory (2006 Launched)

Sik Sik Yuen Biotechnology Mobile Laboratory (BML) Program (2009 Launched)

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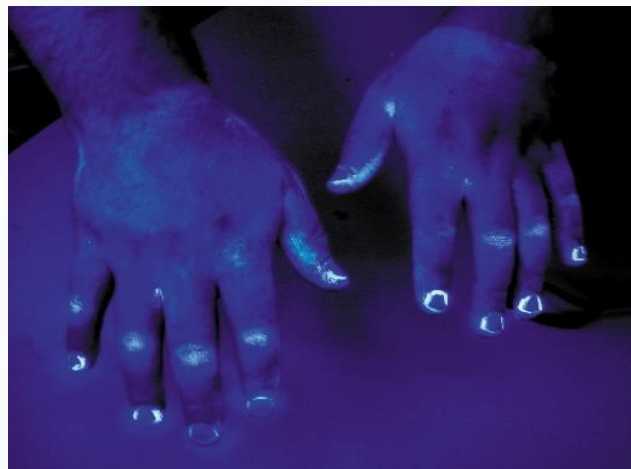
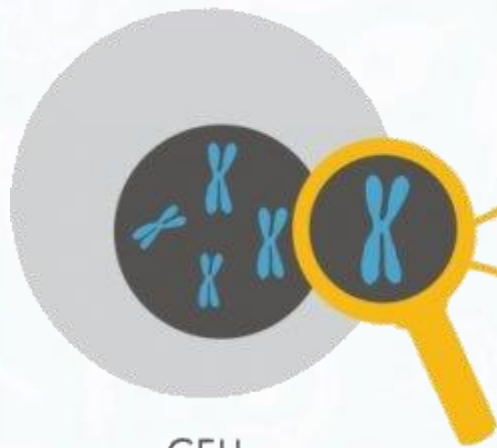
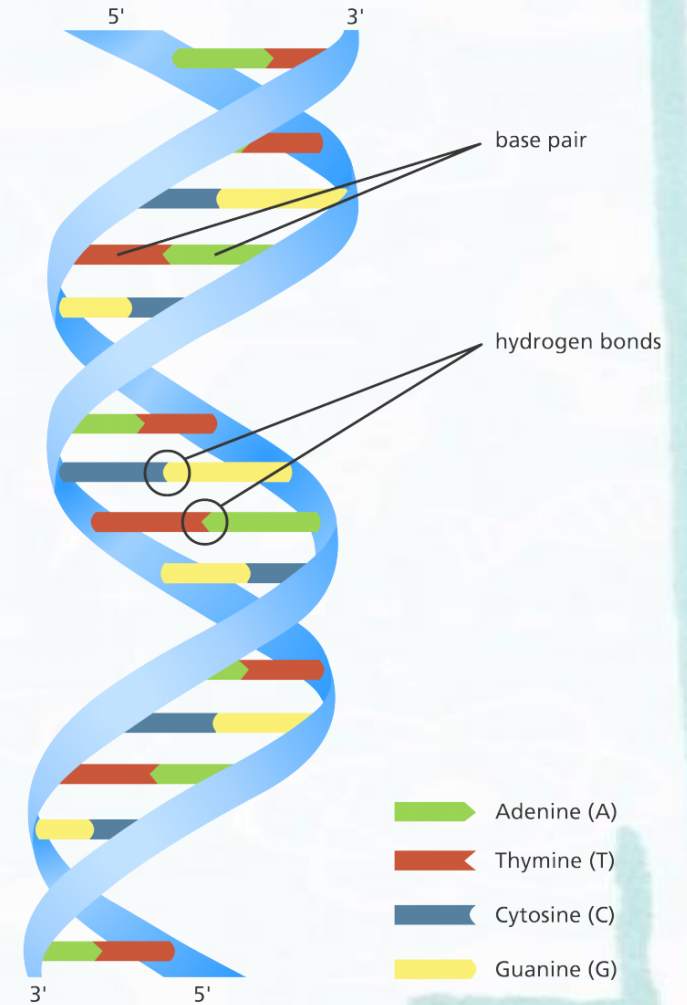
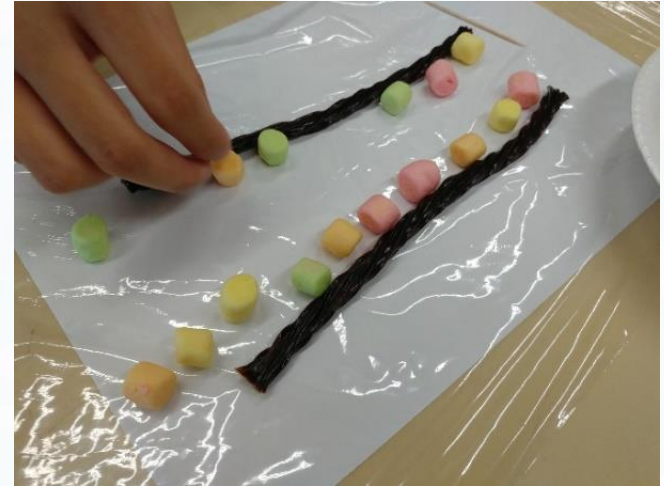
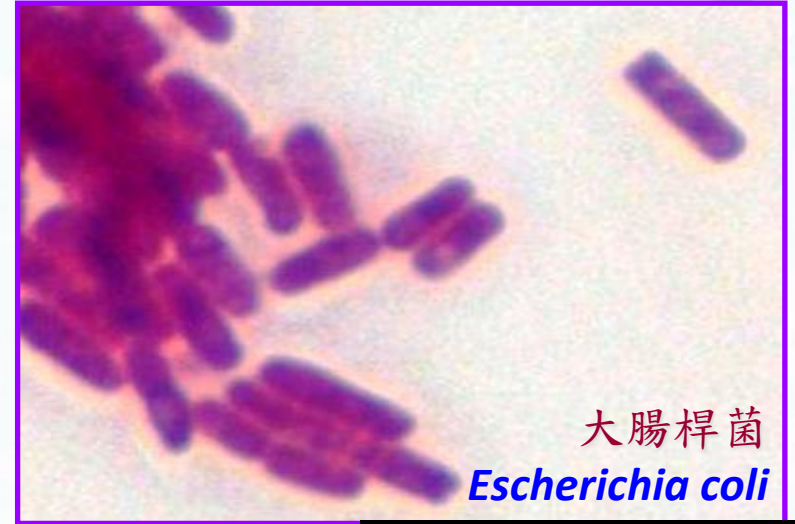
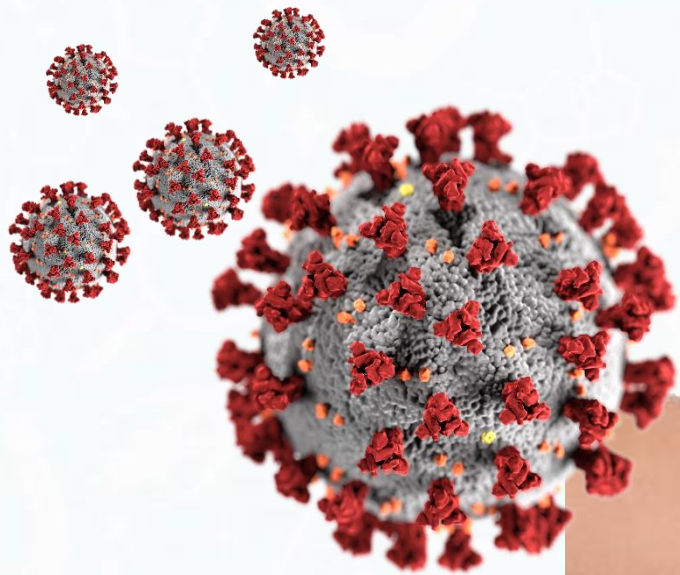
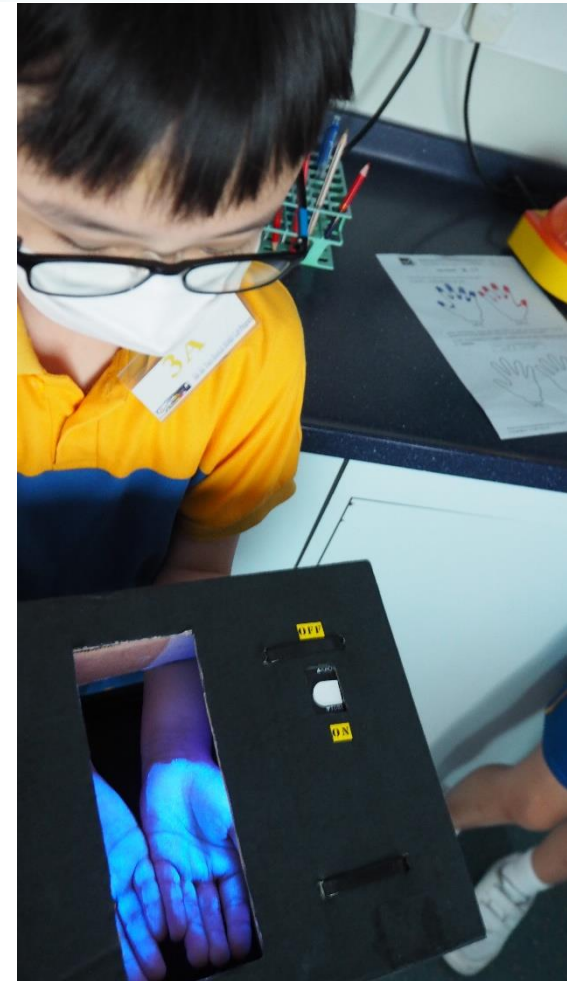
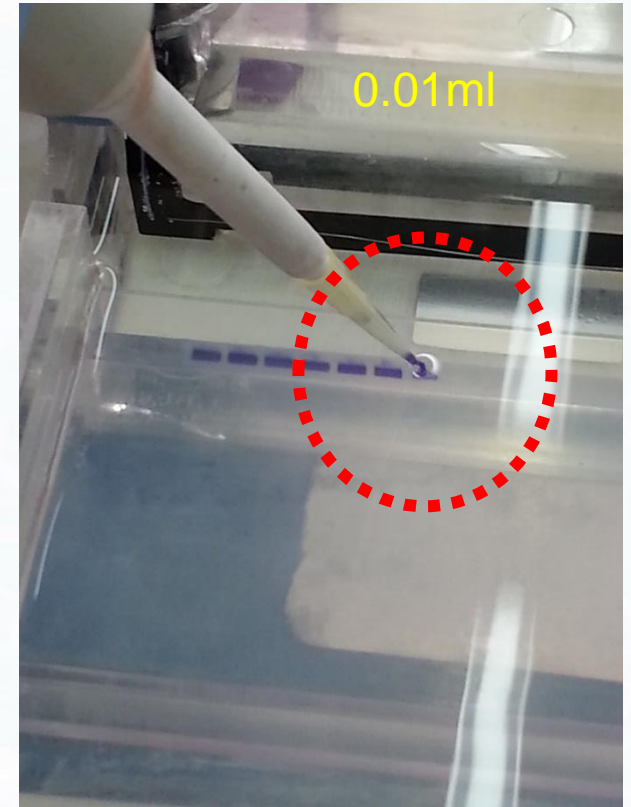
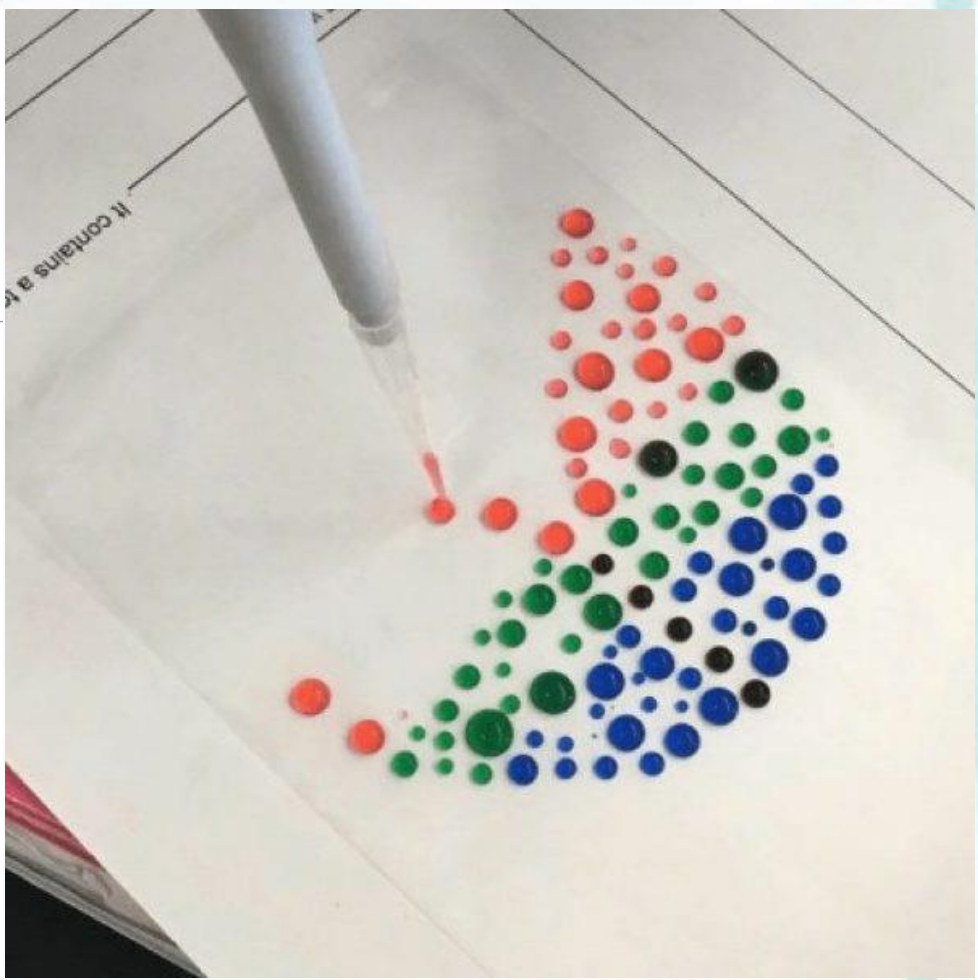
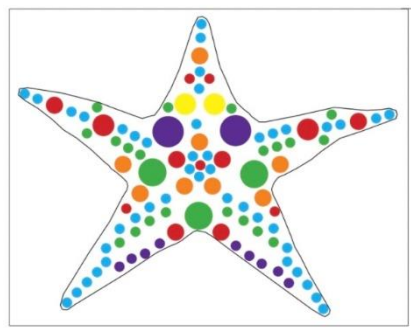
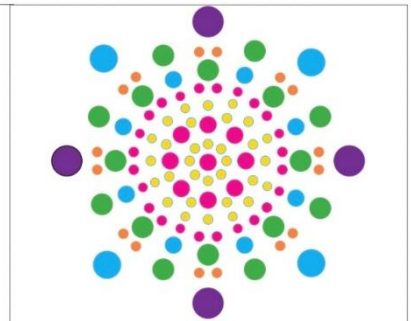
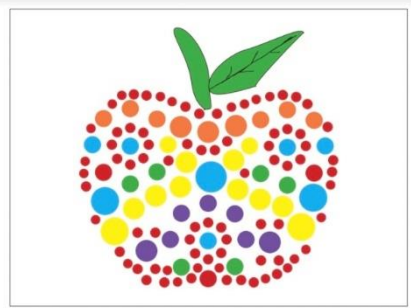
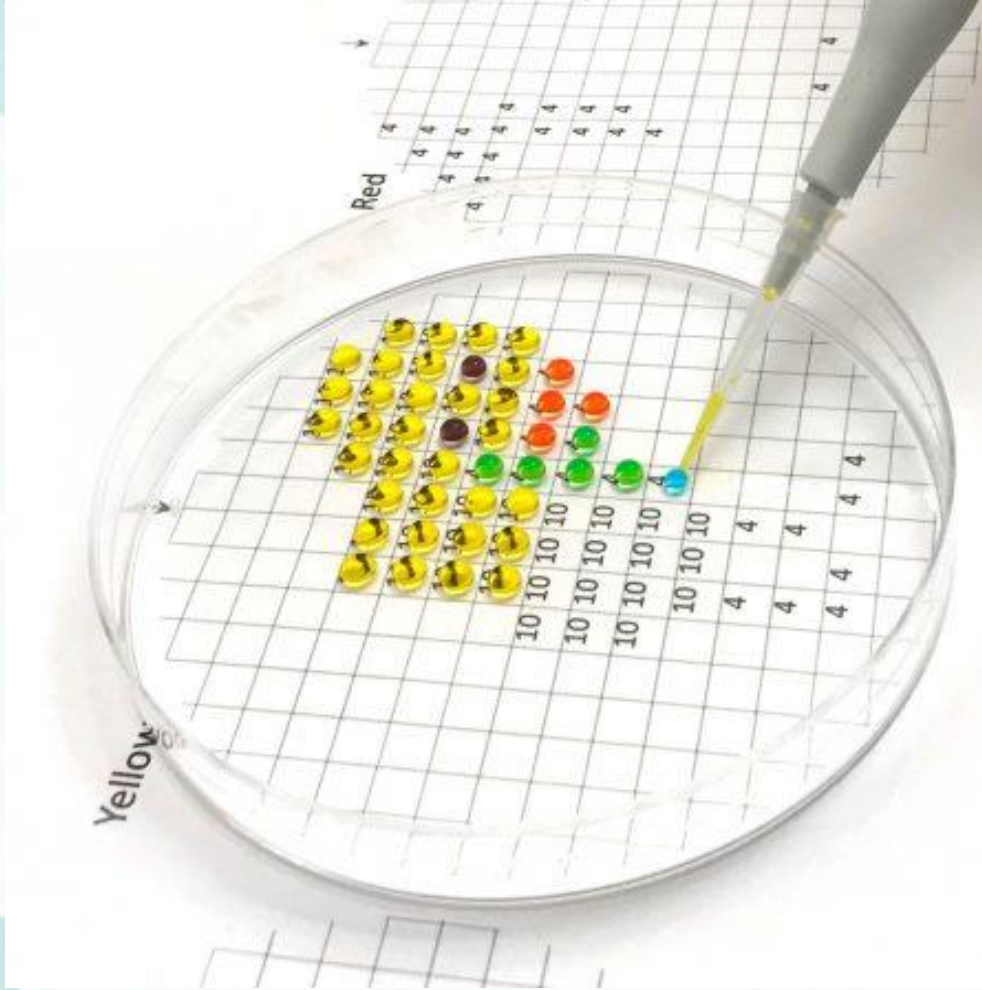
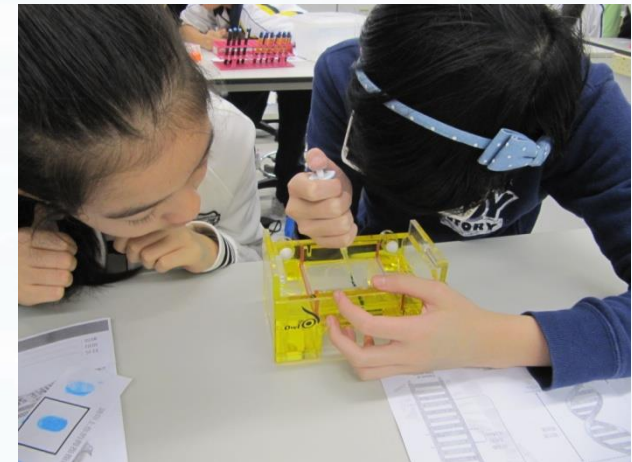
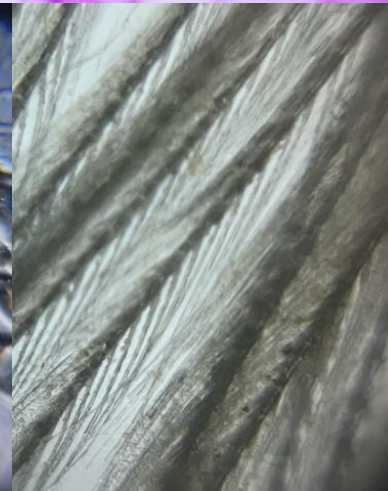
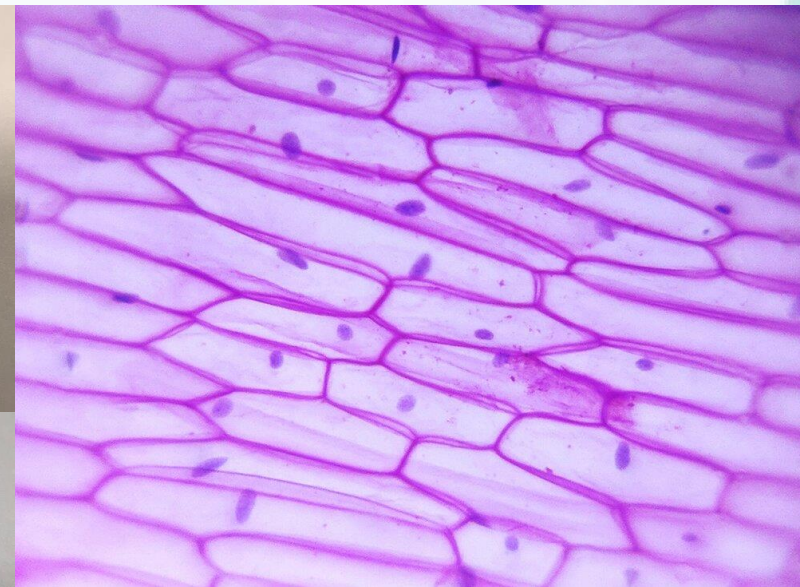
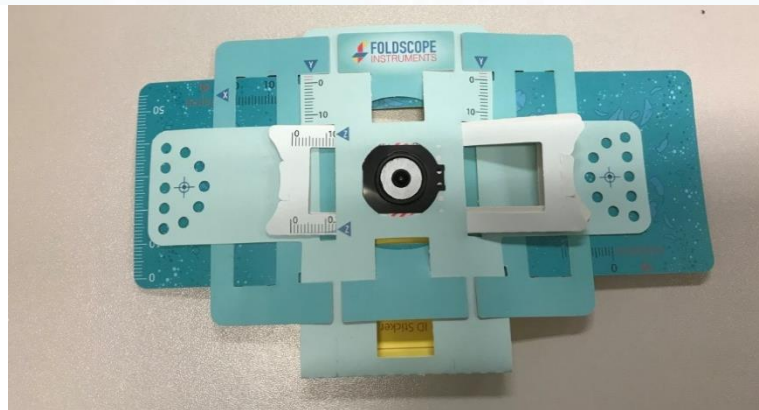
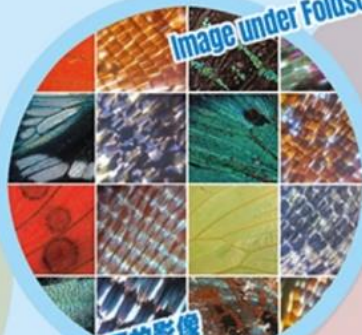
Hong Kong Student Forensic Science Competition 2021

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精彩花絮及比賽結果 Event Highlights & Competition Results

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


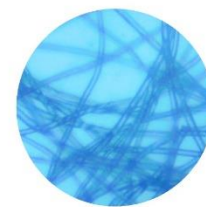
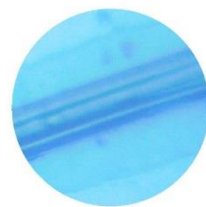
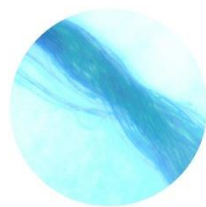
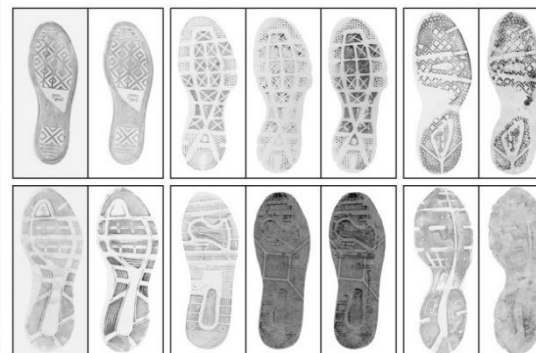
疑犯4
Suspect 4

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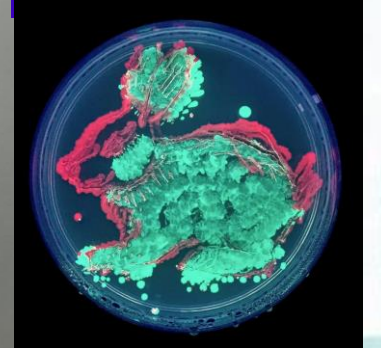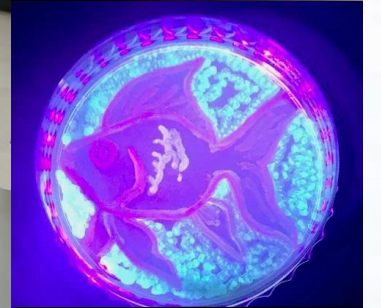
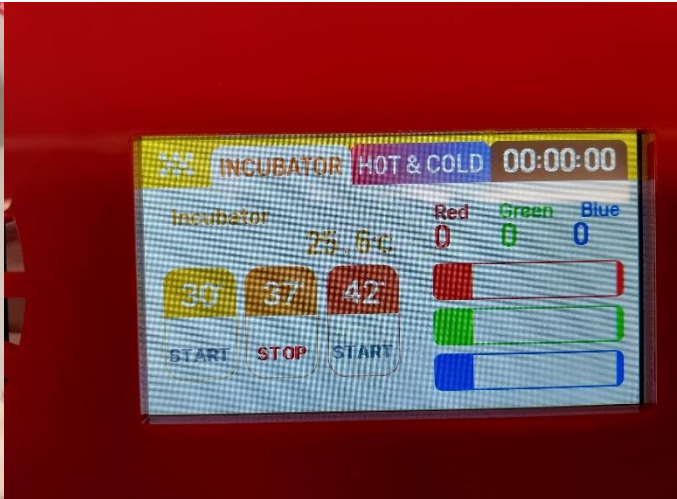
生物工程：轉化 (Transformation)

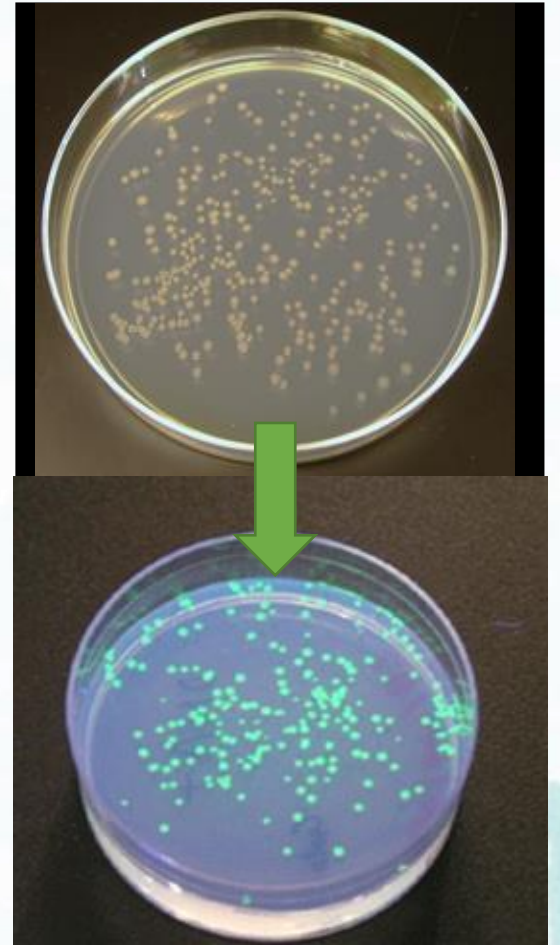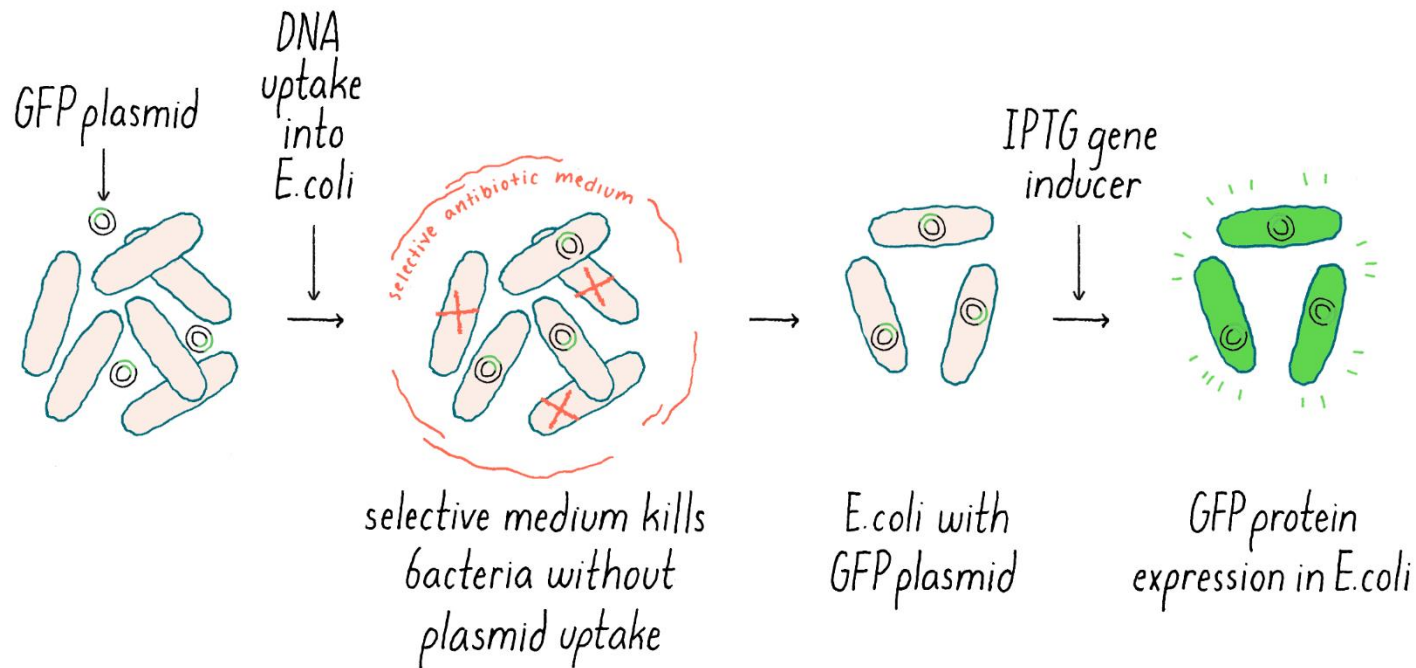TRANSFORMATION AND EXPRESSION OF GFP IN BACTERIA

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
Pre-labs

Pre-lab(s)

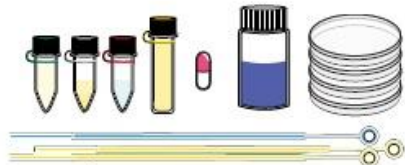Use the Virtual Bioengineer to prepare for the hands-on experience with theory and procedures (30 minutes)

Engineer-it Kit

Day 1

Create petri dishes, LB agar, streak blank bacteria and incubate (45 minutes)

Day 2

Make cells chemically competent & transform with DNA plasmid. Start recovery (45 minutes)

Day 3

Plate transformed cells and incubate (45 minutes)




Day 4

View colonies and analyze results. (10-45 minutes)


Incubation
12-24 hours

Recovery
12-24 hours

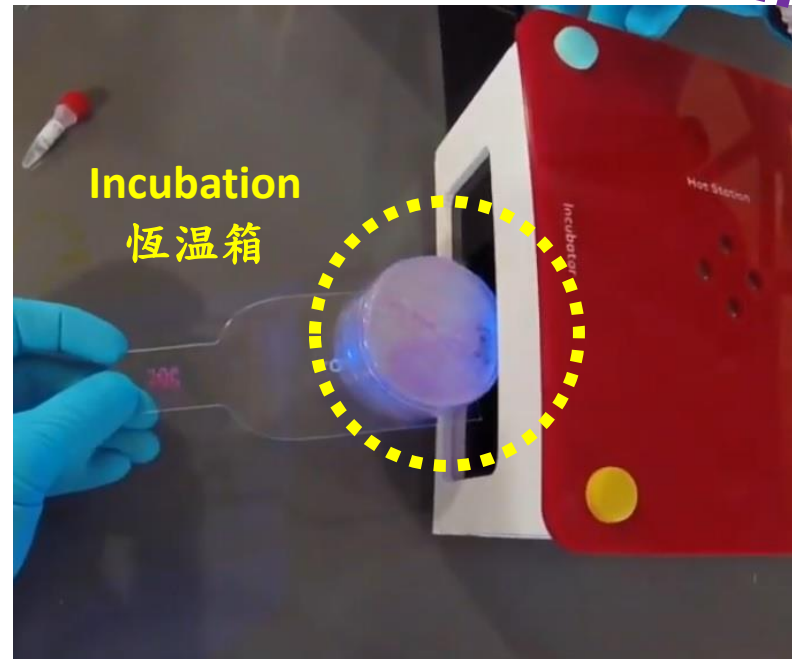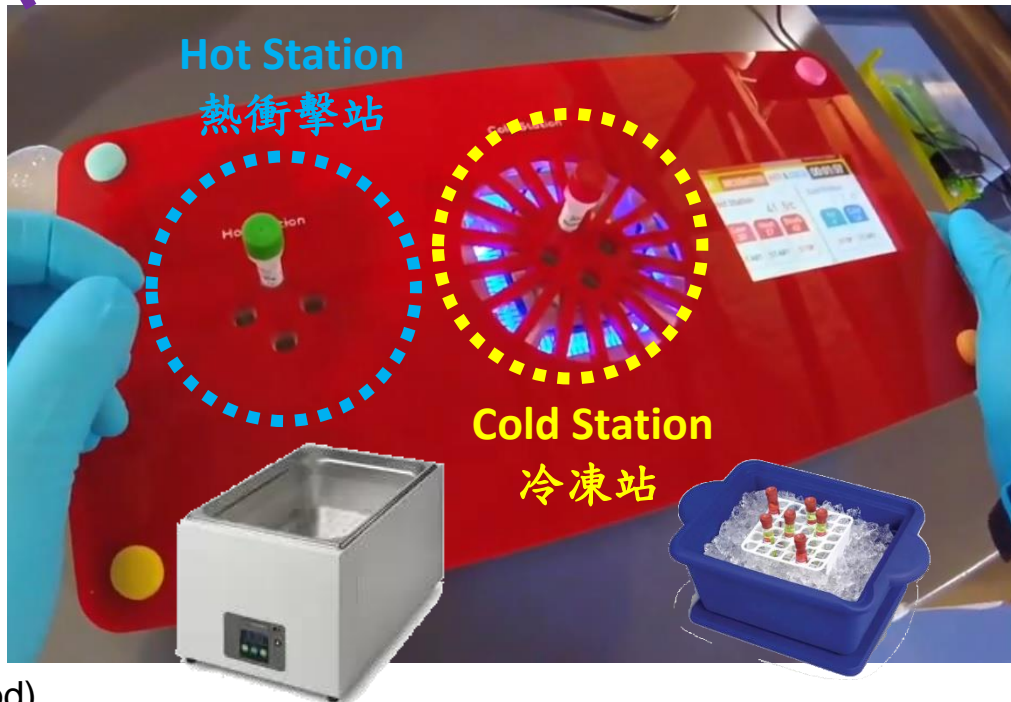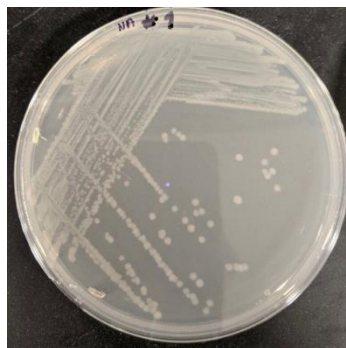Incubation
24-48 hours

劃線法 (streak-plate method)

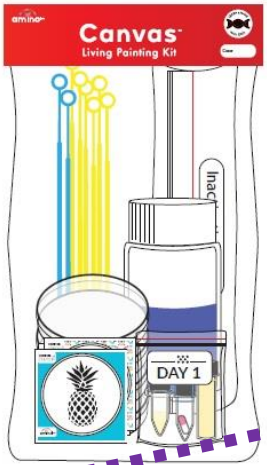Canvas Kit

Day 1

Create painting palette(s)
 Each student group or individual makes LB agar plates with antibiotics (selective plates) and streaks each of the colored bacteria paints from the tubes included in the kit onto one of these selective plates. This amplifies the colorful bacteria so there is enough to paint with. Incubate it so that it becomes the painting palette. **(60 minutes)**


Day 2

Paint your bioart
 Each student or individual can use a blank stencil to draw their picture that will become the bioart. Set the stencil under the petri dish in order to trace it using the bacteria paintbrushes and the painting palette. Incubate the petri dish(es). **(30+ minutes)**

Day 3

See your bioart
 View the living art grow and change color over the next 24-72 hours. Use natural and UV light to see the different colors and photographs! **(10+ minutes)**

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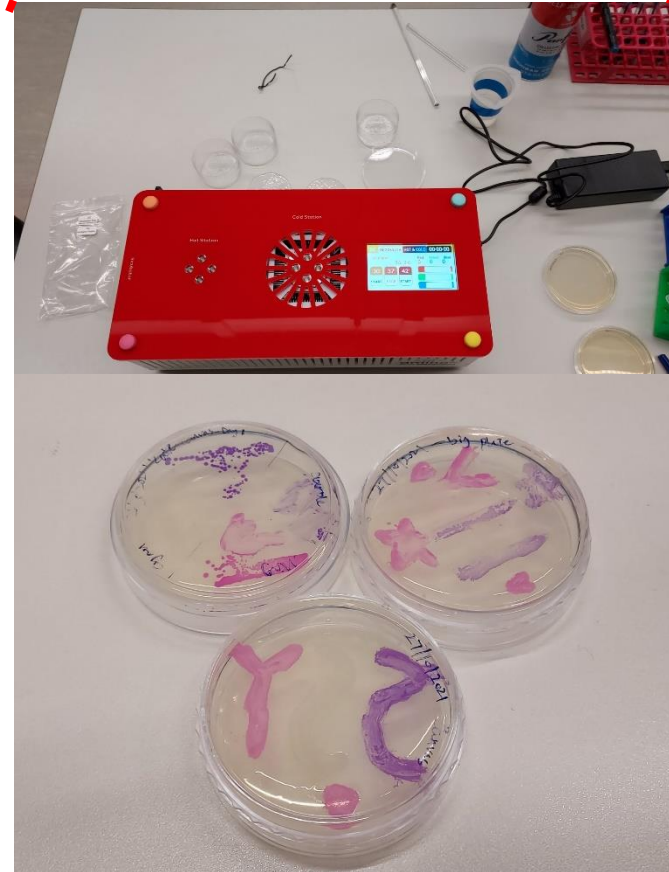Living art by Flowery Branch High School

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This video showcases the QX600™ Droplet Digital™ PCR System which enables advanced six-color multiplexing, allowing clear discrimination of multiple targets. The QX600 System is designed for researchers who need to quantify multiple targets with high accuracy, reproducibility, and sensitivity.

Learn more about QX600™ Droplet Digital™ PCR System at ...

Digital PCR ▶ 全部播放

Bio-Rad's Droplet Digital PCR™ (ddPCR™) systems provide absolute quantification of target DNA or RNA molecules with EvaGreen or probe-based digital PCR assays without using a standard curve. ...

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Edvotek was founded in 1987 as the first company focused on translating cutting-edge biotechnology for the teaching classroom. We work with educators all over the world to demystify science and foster the next generation of scientists through hands-on, active learning activities.

Our equipment is tough enough to stand up to daily use in the ...

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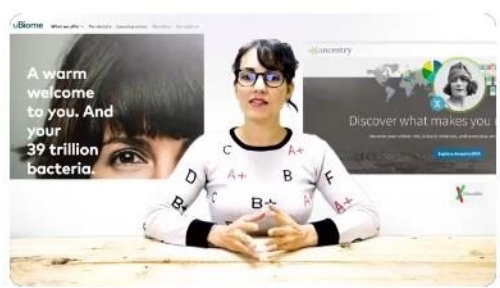
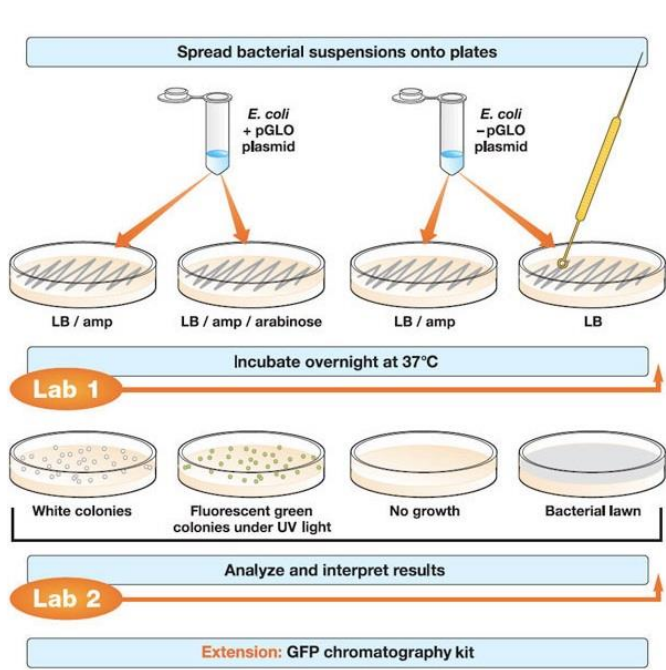
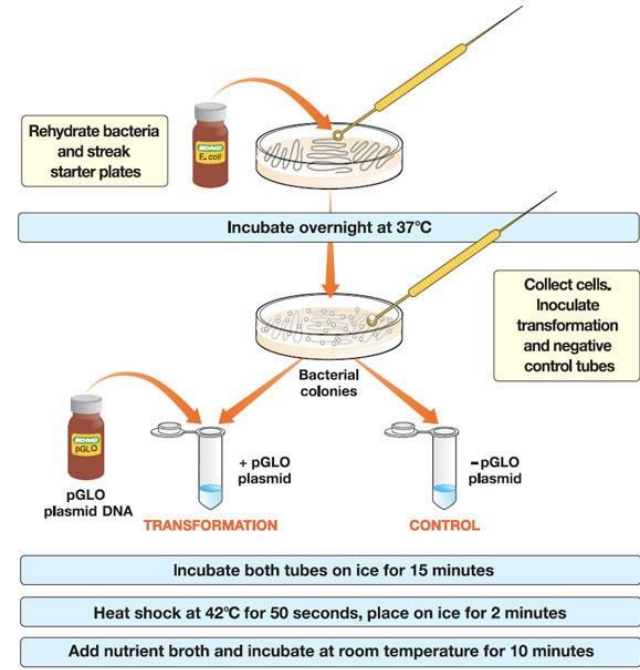
Introducing Foldscope: The Paper Microscope
 觀看次數：8,117次 · 2年前

Foldscope magnifies curiosity - the sense of wonder of the world all around us. From your kitchen sink to the water in the creek running in the backyard; from life forms bizarre and wild to the smallest bacteria crawling in your kitchen sink. See and experience science for yourself! All microscopy clips in the video are taken with a Foldscope.

...

Welcome to the Biorevolution

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Biology-as-a-Technology (BAAT) is the mindset where biology is not just a natural force, but rather a powerful technology platform. As genetic engineering and BAAT will become widespread over the coming decades, there is not better time to start learning about it than now. Become "bio-literate".

Tutorials ▶ 全部播放

- EXTRACTING DNA FROM A STRAWBERRY (6:54)
- Canvas Kit for living (1:16:44)
- The Engineer-it Kit for (1:40:49)
- Smell-it Kit experiment video: (39:34)
- How to make a DIY incubator (12:5...)

Virtual Bioengineer - Canvas Kit Edition

Remember that for any of the Virtual Bioengineer simulators, **you need to be using a desktop computer** as they are not optimized for touch devices like tablets and phones. Go full screen for the best experience!

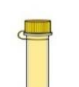
Canvas Kit

1. Make LB Agar Plates
2. Grow Paint Palette
3. Incubate
4. Paint Living Canvas
5. Incubate

Materials

Sterile Water

LB Agar Powder

Microwave

Antibiotics

Nitrile Gloves

Petri dishes

Start Making LB Agar Plates

YouTube

DAY 1 ADD AGAR

2:13 / 13:00

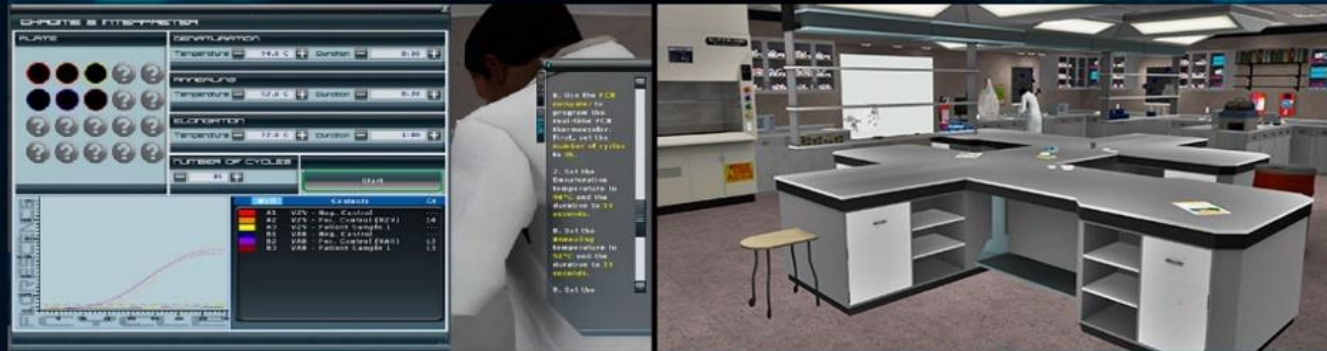
Zero to Genetic Engineering Hero: Create Living Paintings with Canvas Kits

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MBT Mission Biotech

Free download. Hours and hours of engaging, immersive, 3D gameplay.

In response to the COVID-19 outbreak, we are making Mission Biotech FREE to the public. This fun, immersive game contains real-world laboratory work in the field of biotechnology, targeted for middle schoolers and above. Includes techniques used by real-world medical workers to test for the currently circulating coronavirus including DNA Extraction and PCR. Funded by NSF, this game encourages everyone to explore careers in biotechnology and work against the clock to identify the threat!

For PC/Windows only

PLAY FOR
FREE

UF UNIVERSITY of FLORIDA

Number of downloads: 1700

Download [Reference Materials](#)

Download [Gameplay Images](#)

Mission Biotech

User Guide


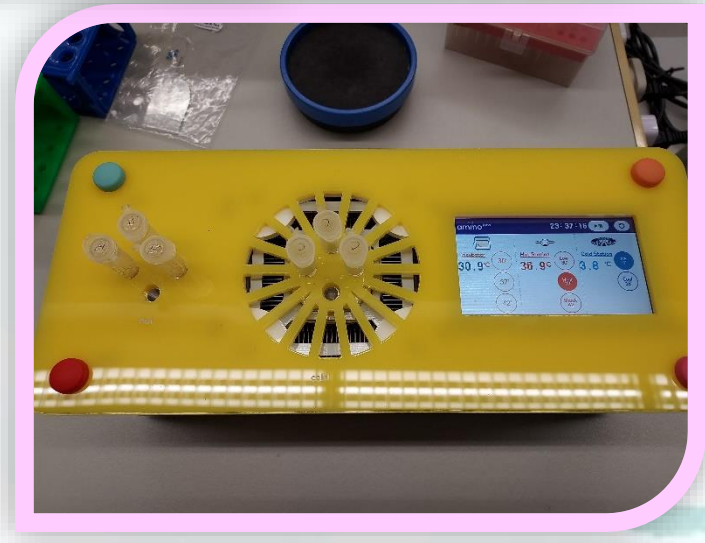
VIRTUALHEROES
A DIVISION OF APPLIED RESEARCH ASSOCIATES, INC.



Table of Contents

| | |
|-------------------------------------------|-----------|
| 1. Overview | 4 |
| 1.1 Goals | 4 |
| 1.2 Mode of Play | 4 |
| 1.3 Features | 4 |
| 1.4 Platform | 4 |
| 1.5 Minimum System Requirements | 4 |
| 1.6 Recommended System Requirements | 5 |
| 2. Objective | 5 |
| 3. Launching the Game | 5 |
| 4. Mission Details | 5 |
| 4.1 Mission Progression | 5 |
| 4.2 Adventure Interactions | 6 |
| 4.3 Unlocks | 6 |
| 4.4 Tests | 6 |
| 5. Movement and Control | 6 |
| 6. Character Selection | 7 |
| 7. Badge Creation | 9 |
| 7.1 Name Character | 9 |
| 7.2 Choose Badge Style | 9 |
| 7.2.1 Rank | 9 |
| 7.2.2 Achievements | 9 |
| 8. User Interface | 10 |
| 8.1 User Interface Elements | 10 |
| 8.1.1 Cursor | 10 |
| 8.1.2 Inventory | 11 |
| 8.1.3 Active Item | 11 |
| 8.1.4 Character Badge | 12 |
| 8.1.5 Map | 12 |
| 8.1.6 Protocols Cheat Sheet | 12 |
| 8.2 Software Menus | 12 |
| 8.2.1 Main Screen | 12 |
| 8.2.2 Options Screen | 12 |

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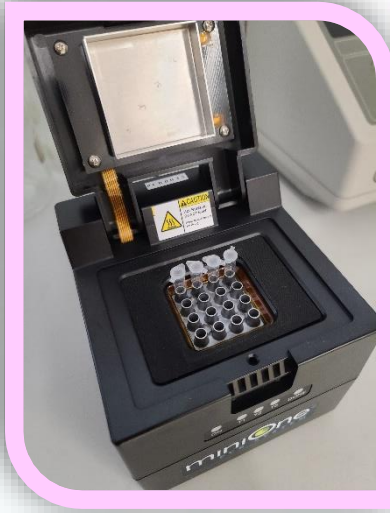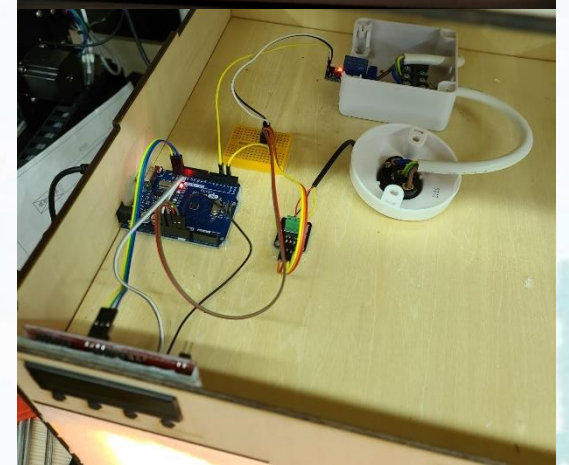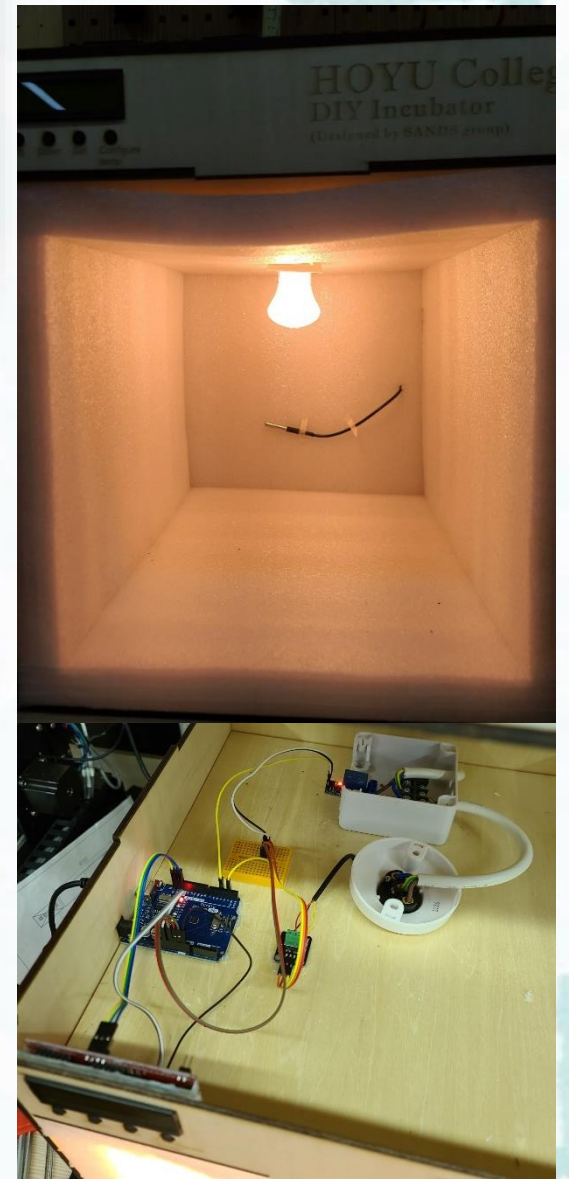SIK SIK YUEN

Biotechnology Mobile Laboratory Program

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